



Migrating to a new PC

On the old PC


- Insert a memory stick into an available port
- Select 'Data Management', then 'General'

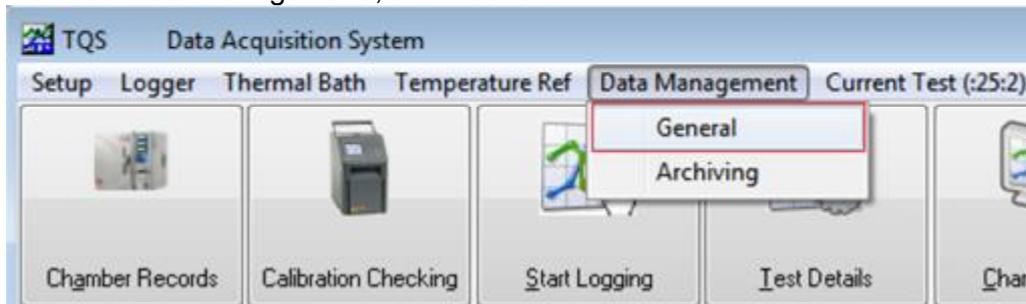

- Select the correct drive for the memory stick on the right side

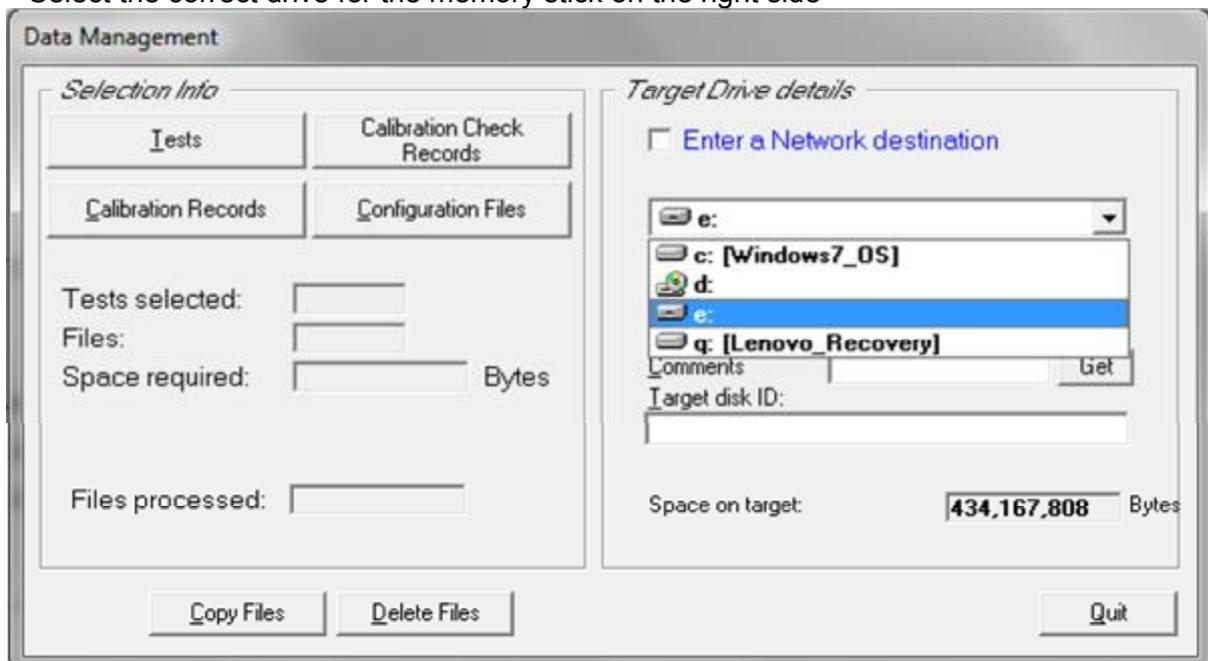

- Select 'Configuration Files'

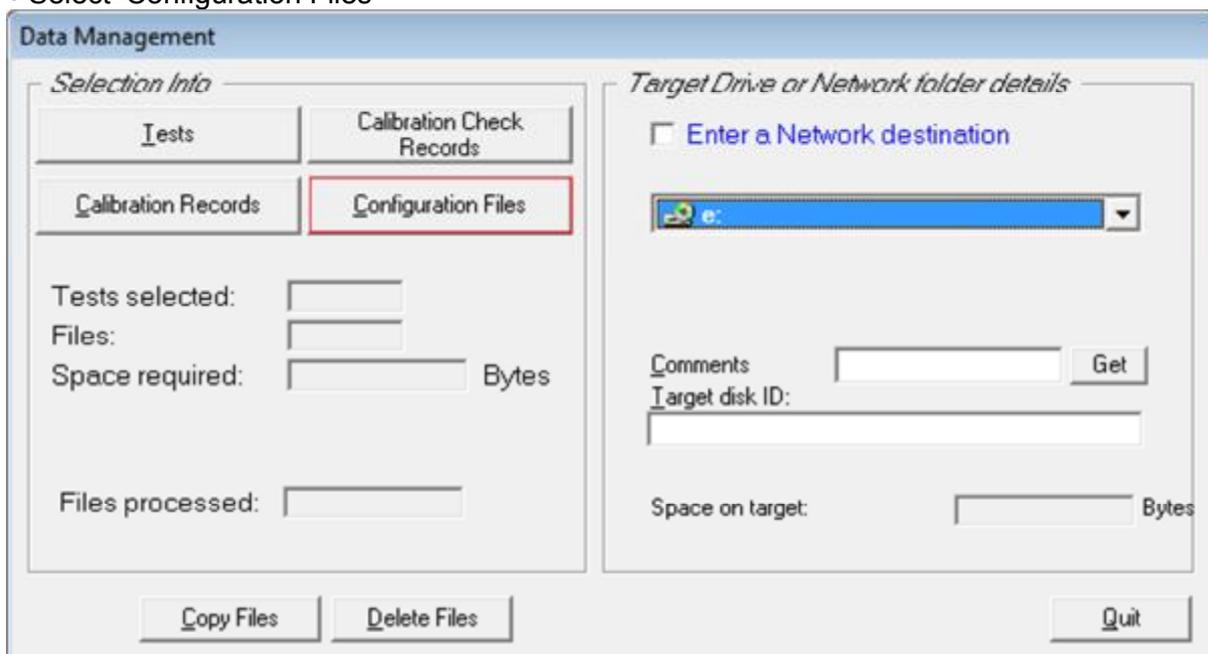

- Choose which files you wish to copy across
Basic Settings – which logger, heat bath is set to default etc


Channel Configurations - including labels etc

Security Settings – user IDs and passwords

Calibration Files – historic calibration data

Chamber Records – copies chamber records (but not historic data)

Test Specifications – including any user defined Test Specs

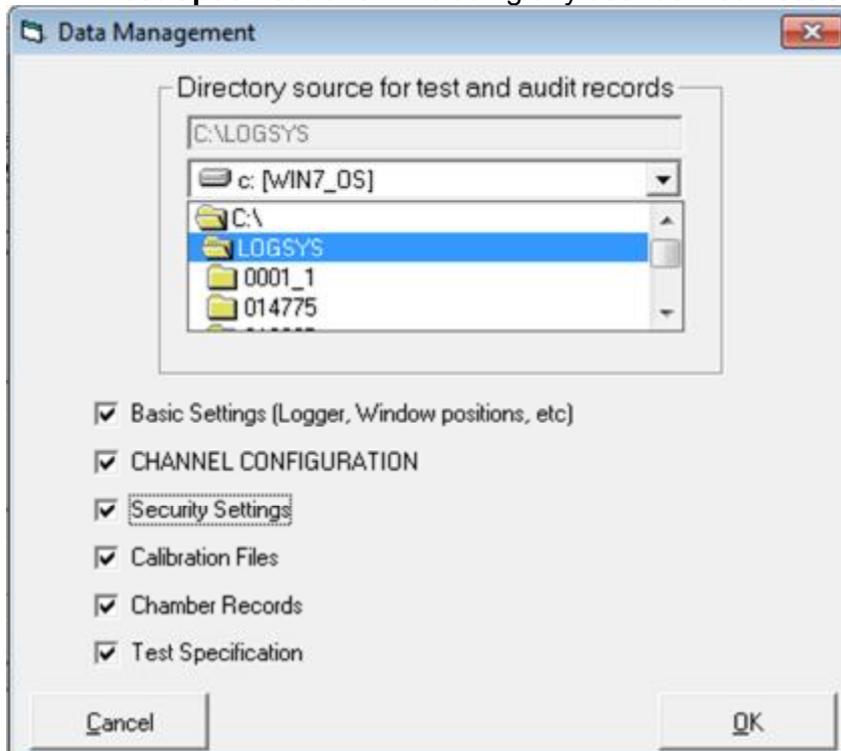

- Press OK. A folder called 'Logsys will be created on your memory stick.
- Safely eject the memory stick from the first PC, and insert it into a spare slot on the destination PC
- Copy the Logsys folder from your memory card and paste into your C:/ drive. When warned that a file with this name already exists, choose 'Copy & Replace'